



Material Characterization of Rock Sample using Thermal Vaporization/Pyrolysis-GC/MS Technique

Authors:

Dr. Tom Malloy & Dr. Adry Bissada, University of Houston, Center for Petroleum Geochemistry
Rojin Belganeh, Frontier Labs Americas

Introduction

One of the first steps in the standard protocol for geochemical characterization of rock samples is Rock-Eval analysis. In its simplest form, this process entails grinding the rock sample and then determining the amount of free or adsorbed matter by heating the sample in an inert gas flow to 300 °C for 3 minutes (S1) and then ramping the temperature to 650 °C @ 25 °C/min (S2) to pyrolyze the heavier material. The hydrocarbons are measured with an FID and CO₂ (S3) by an infrared detector. The value of S1 is related to the recoverable oil, and S2 of the potential for oil production. A measure of the maturity may be derived from the temperature of the maximum in the S2 curve T_{ps2} .

Key Words

- Rock-Eval Analysis
- Geochemical Characterization
- Mining
- Oil Production

Method

In the Evolved Gas Analysis (EGA) mode, the Thermal Vaporization / Multi-Shot Pyrolyzer (EGA/PY-3030D) system from Frontier Laboratories, paired with a Frontier AS-1020E autosampler, a Frontier SS-1010E Selective Sampler, and a Frontier MJT-1035E MicroJet Cryo-Trap may be used to determine the basic parameters listed above. This is accomplished by coupling with an Agilent 7890B GC-FID/5977 MSD system and an Agilent 3180B two-way splitter with makeup gas. In this configuration, an inert Ultra Alloy capillary is used between the GC injector and the splitter.

It is possible to obtain much more detailed molecular information without extensive solvent extraction to obtain extracted organic material (EOM). Using different modes of operation of the multi-shot pyrolyzer with a GC separation column in place instead of a deactivated EGA tube, this is done in two steps.

In this experiment, a rock sample, referred to as sample A, with a depth of 7845.42 feet is used for material characterization. The ground rock sample is heated to 300 °C and held for 20 minutes while the

evolved gases are cryo-trapped at the head of the column with the MicroJet Cryo-Trap. At that point, the sample is automatically removed out of the furnace to an ambient holding position, the oven and the nitrogen cooling turned off and GC/MS and FID chromatograms are obtained. With the GC/MS operated in SIM/SCAN mode and the split to the FID, three chromatograms with identical retention times, are obtained in a single run. This corresponds to the S1 peak. After this is complete, the nitrogen cooling is turned back on, the sample reintroduced into the oven and the evolved gases are cryo-trapped as before as the oven is heated from 300 °C to 650 °C corresponding to S2. Again, 3 chromatograms are obtained in a single run.

The GC was operated at a constant flow of 1.2 mL min⁻¹ with a temperature program of 70 °C for 3 mins; a ramp of 5 °C per min to 200 °C; a ramp of 4° per min to 320 °C and held at that point for 20 min. The flow was split between the FID and the MSD. The MSD was operated in SIM-SCAN mode.



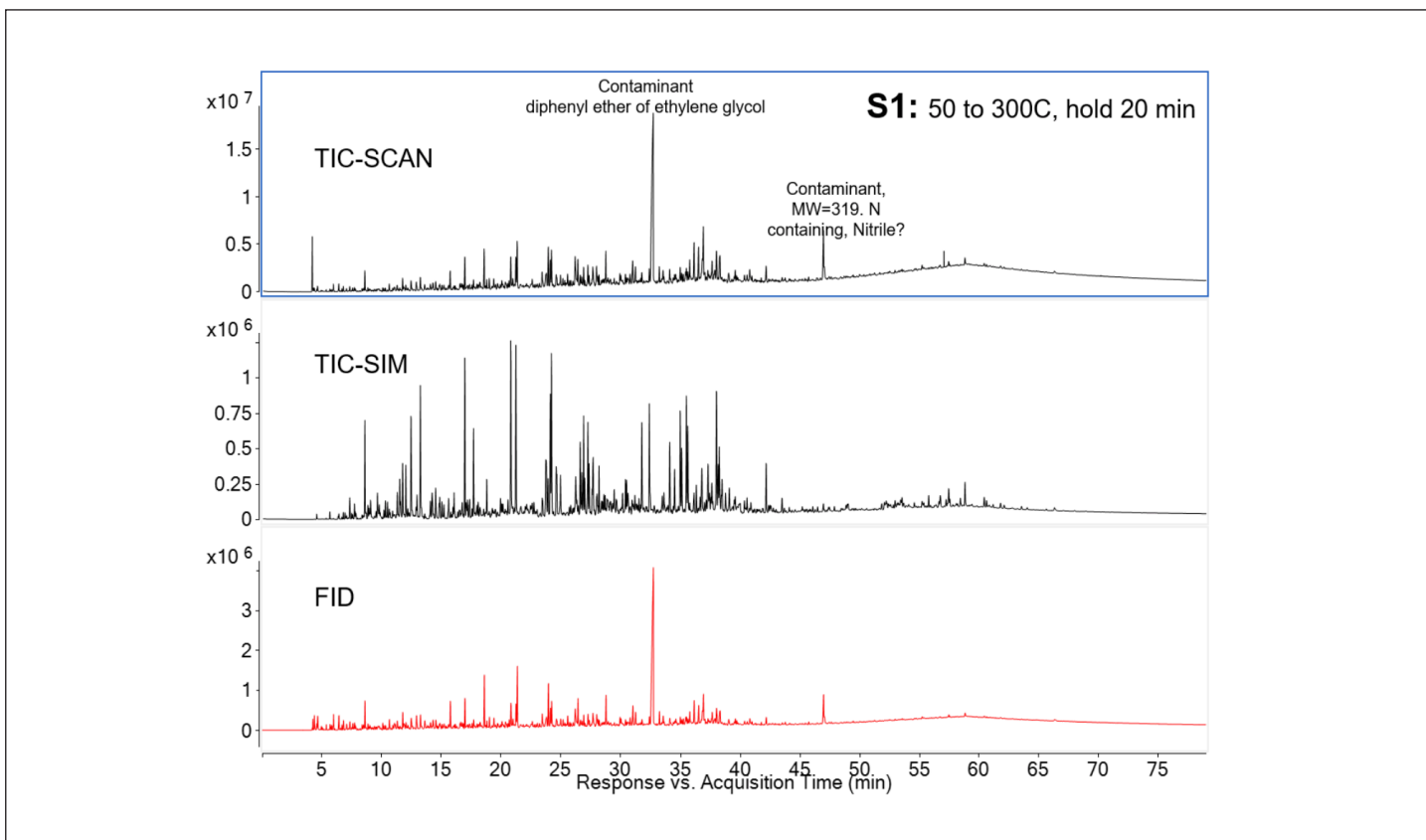


Figure 1. Three Chromatograms Obtained from GC/MS in SIM/SCAN Mode & Split to FID

Results

Figure 1 shows the GC/MS total ion current (TIC) SCAN from $m/z=20$ to 650 (top), the TIC selected ion monitoring (SIM) chromatogram (middle) and the flame ionization detector chromatogram (bottom) all obtained in a single run. Immediately obvious from the TIC SCAN is the presence of two contaminants. The most prominent is the diphenyl ether of ethylene glycol and the second is (actually 2 peaks) nitrogen-containing compound(s) with nominal molecular weight 319. Neither of these will interfere with further analyses.

The Frontier EGA/PY-3030D Multi-Shot Pyrolyzer can analyze any organic materials using some or all of the following techniques:

- Evolved Gas Analysis (EGA)
- Thermal Desorption (TD)
- Heart-Cutting (HC)
- Pyrolysis
- Reactive Pyrolysis (RxPy)
- Single, Double or Multi-Shot Analysis

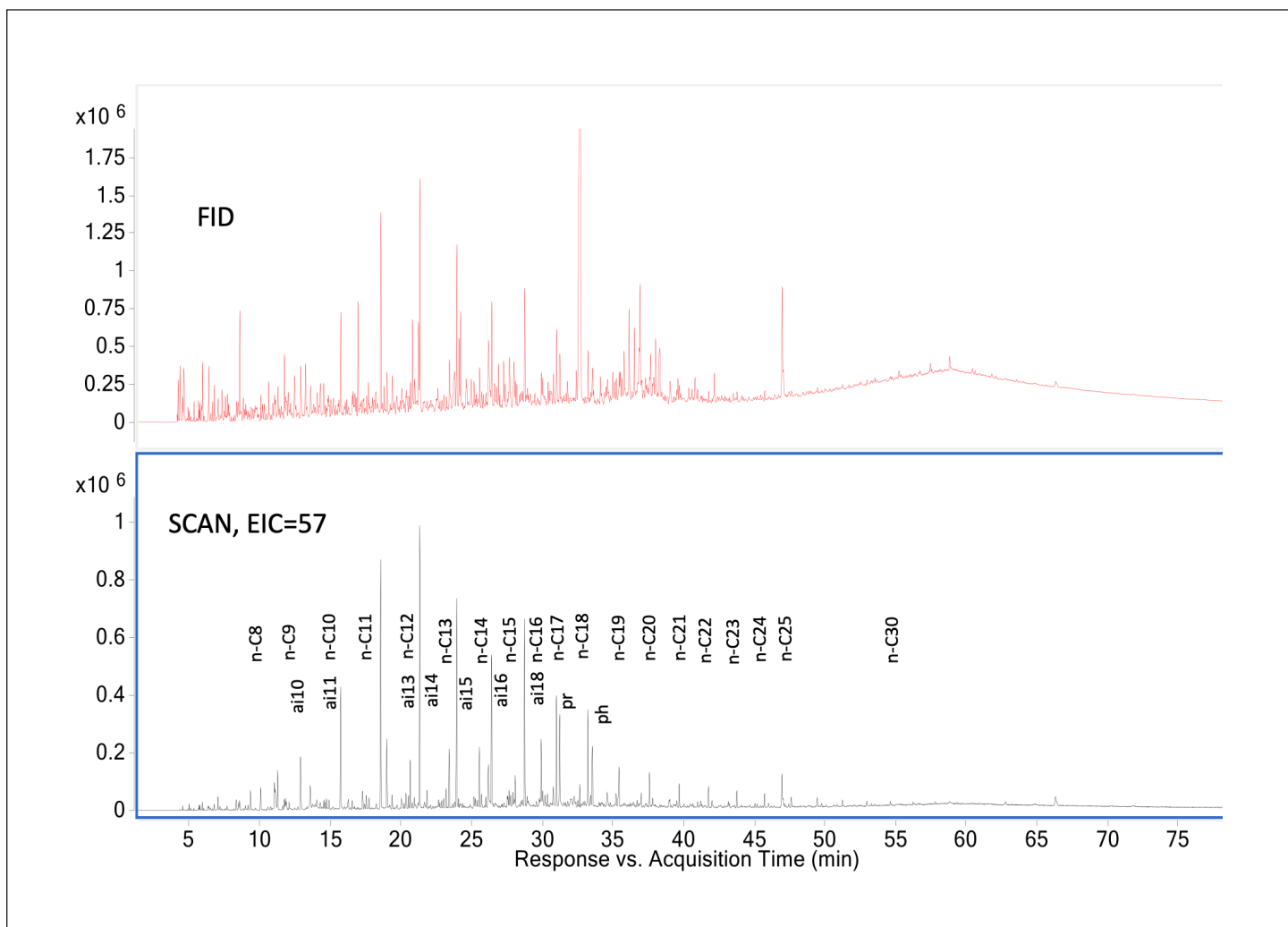


Figure 2. FID & EIC ($m/z = 57$) for n-alkanes and acyclic isoprenoids (ai).

Figure 2 shows the FID trace (top) and an extracted ion chromatogram (EIC) at $m/z = 57$, corresponding primarily to alkanes. Assignments for the n-alkanes and some of the acyclic isoprenoids are indicated in the Figure.

Measurements of peak areas from the FID chromatogram, along with an expanded plot of the FID trace from 4 to 9 minutes are given in **Table 1** in **Appendix A**.

Figures 3 and 4 are EIC plots for substituted benzenes, naphthalenes, phenanthrenes, and dibenzothiophenes. The distribution of isomers and ratios among the alkanes and acyclic

isoprenoids as well as the various aromatics are used to derive geochemical properties, including thermal maturity. Measurements of SIM areas/heights are used for the aromatics.

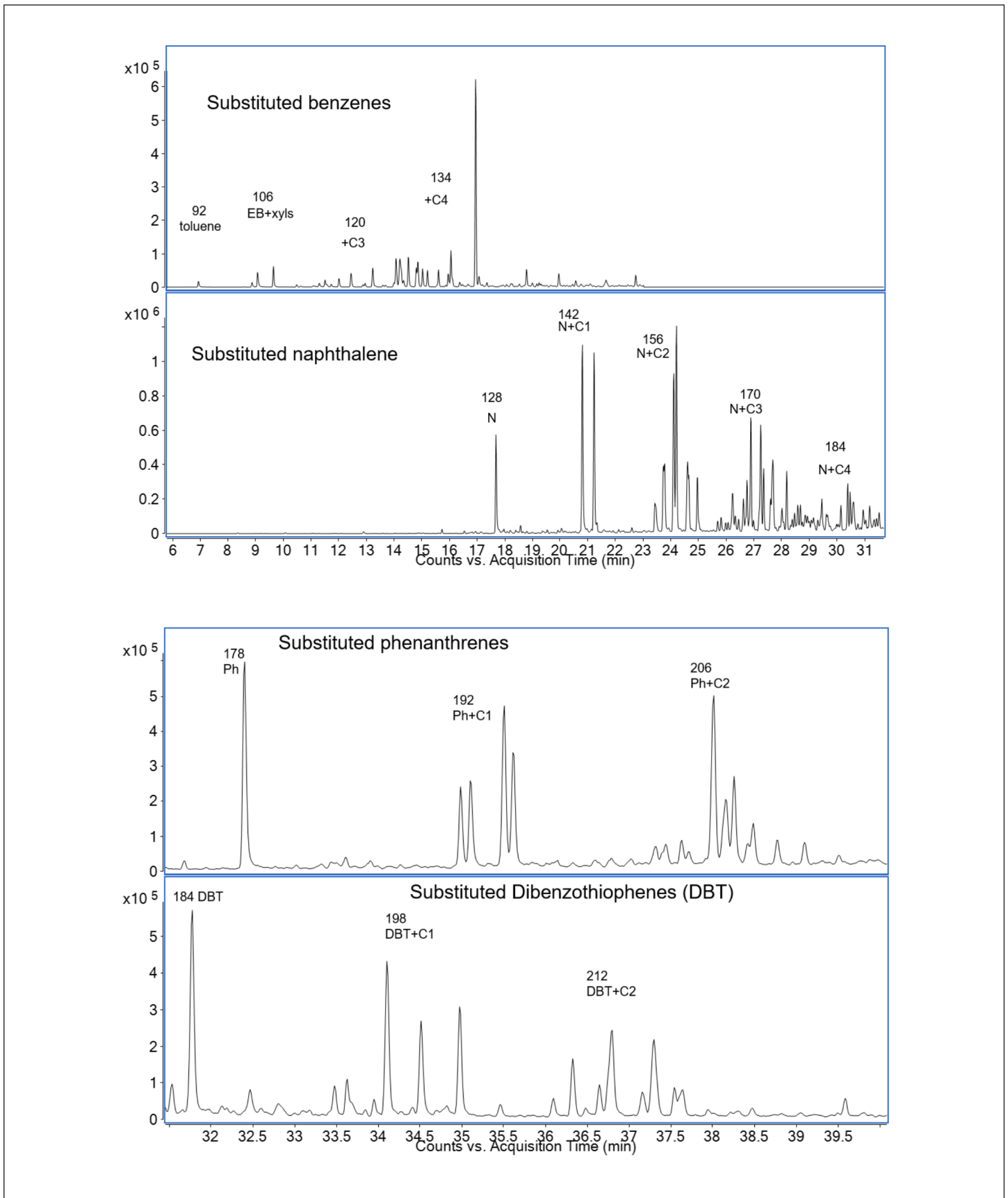


Figure 3 & 4. EIC for Substituted Benzenes, Naphthalene, Phenanthrene, and Dibenzothiophenes.

Even more diagnostic information is obtained from the measurement of biomarkers; terpanes and hopanes, steranes, and mono- and di-aromatic

steroids. **Figure 5** shows the fragmentograms for $m/z = 191.2$, 217.2 and 218.2 . The first corresponds to terpanes and hopanes and the last two to steranes.

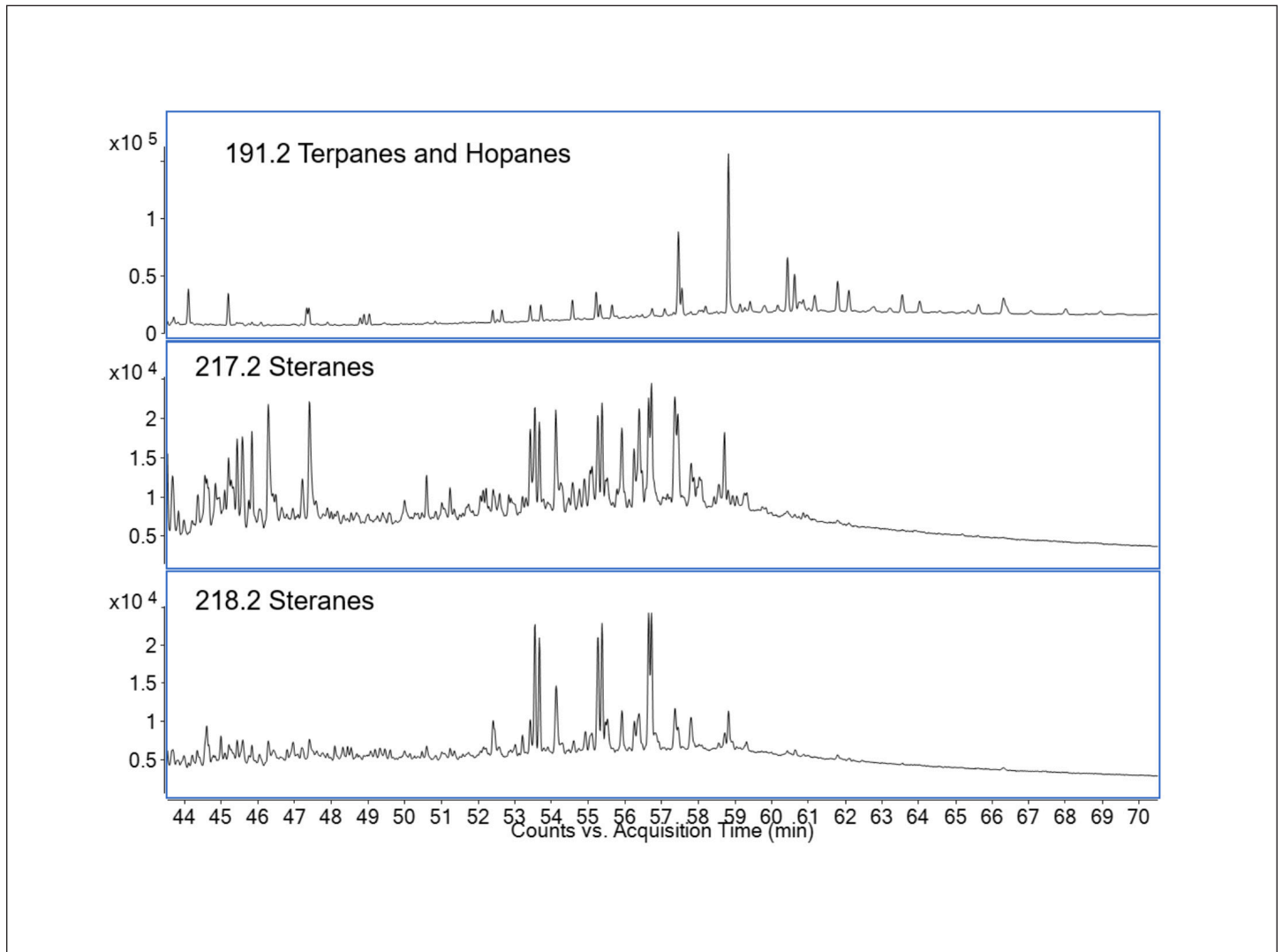


Figure 5. Extracted Ion Chromatograms for Biomarkers (Terpanes, Hopanes, and Steranes).

Figure 6 below shows the fragmentograms for $m/z = 231.1$ and 253.2 , the tri-aromatic and mono-aromatic steroids.

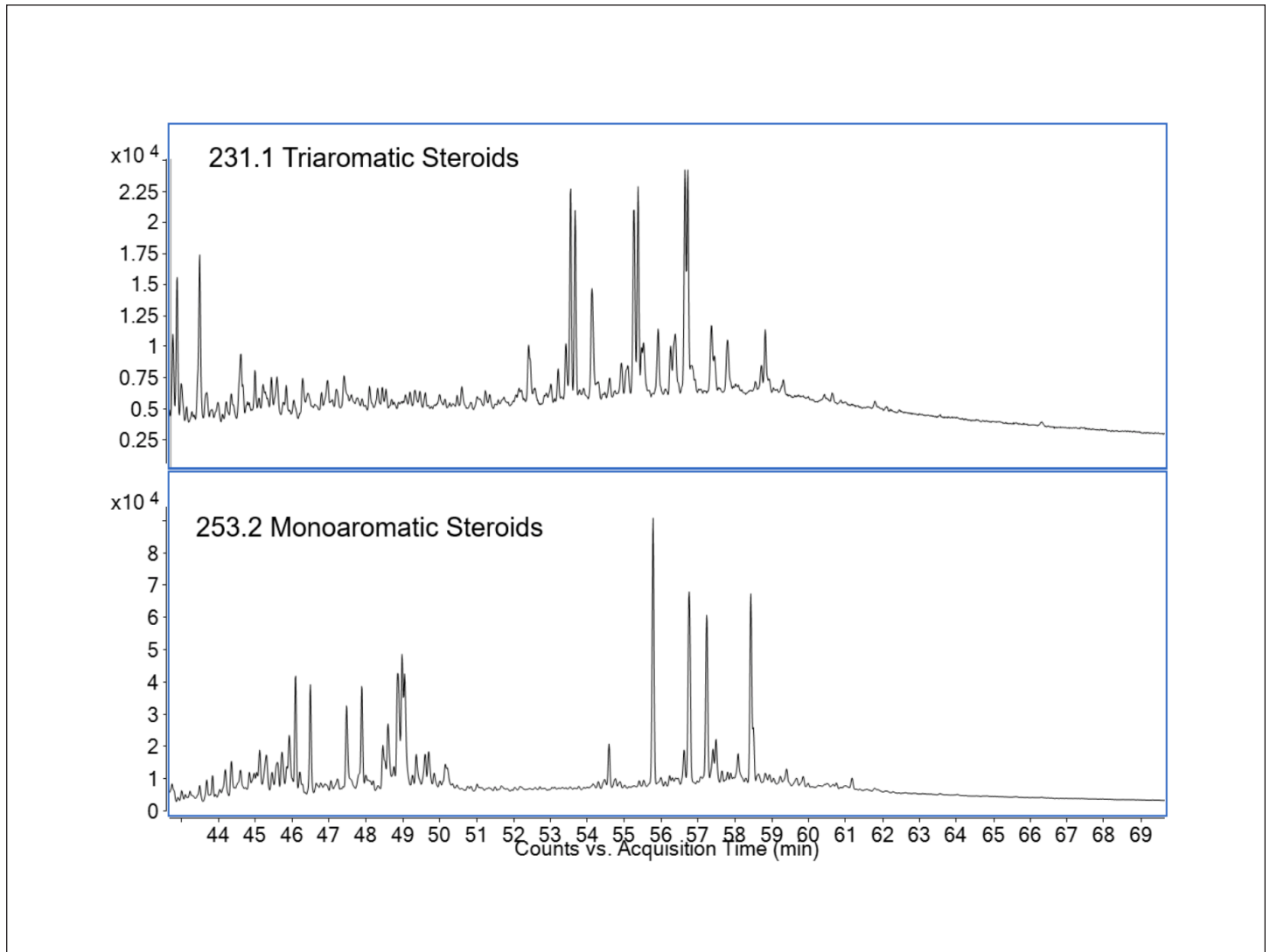


Figure 6. Tri-aromatic & Mono-aromatic Steroids.

Table 2 shows the results for the terpanes and hopanes relative to 30-Hopane normalized to 100 (**Appendix B**). The presence or absence of specific biomarkers and ratios of combinations of biomarkers

yields information on the conditions during deposition, age, maturity, and sometimes lithology.

Figure 7 below is a different sample with assignments shown on the Figure for illustration.

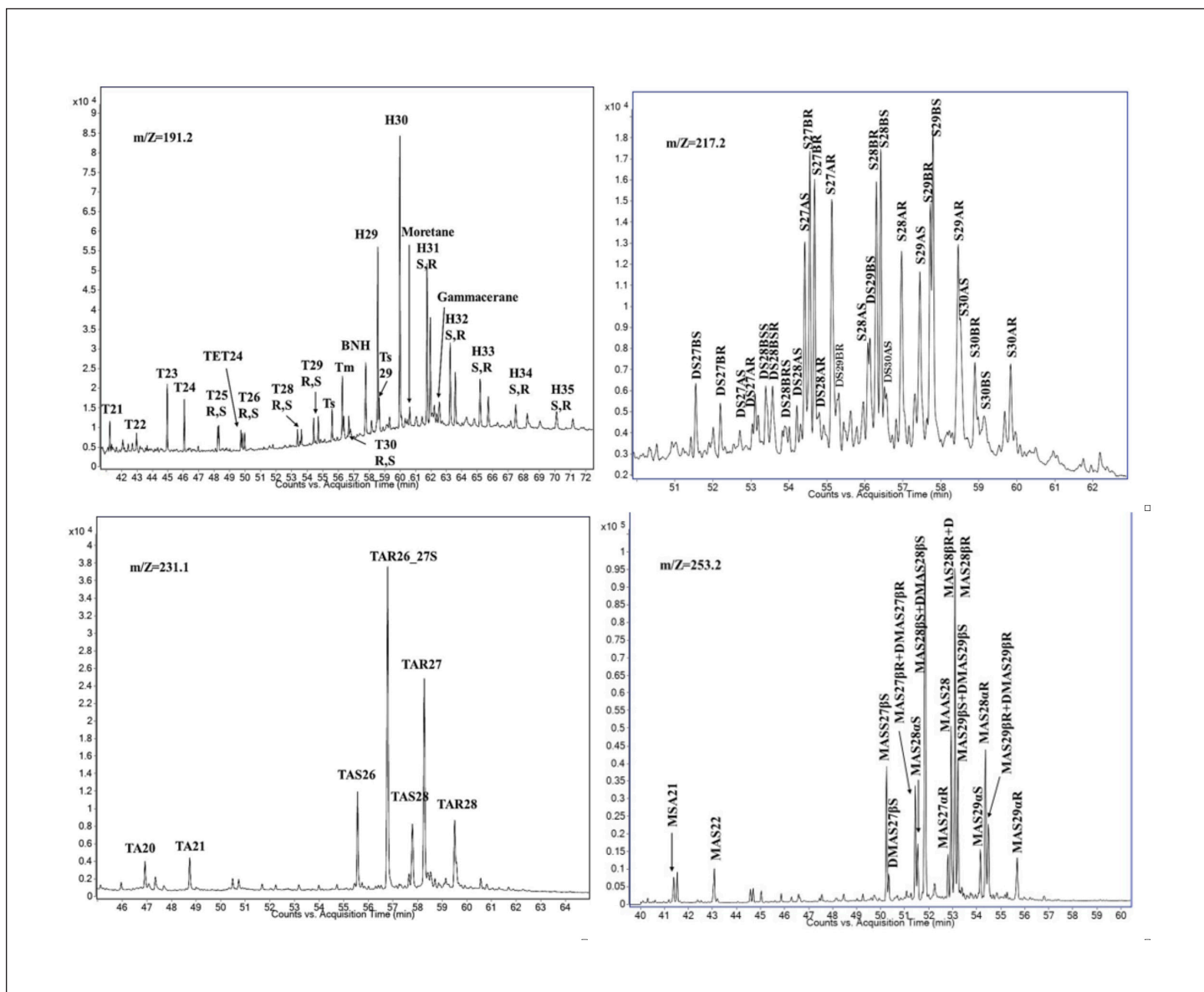


Figure 7. Illustrating Figure for Peak Assignments.

Figure 8 shows the three chromatograms collected for the S2 peak: the TIC SCAN, the TIC SIM and the FID. The TIC SCAN is used along with a library to identify

peaks that may be in question. Many small alkenes are observed early on. The progression of alkane peaks is accompanied by the corresponding alkene.

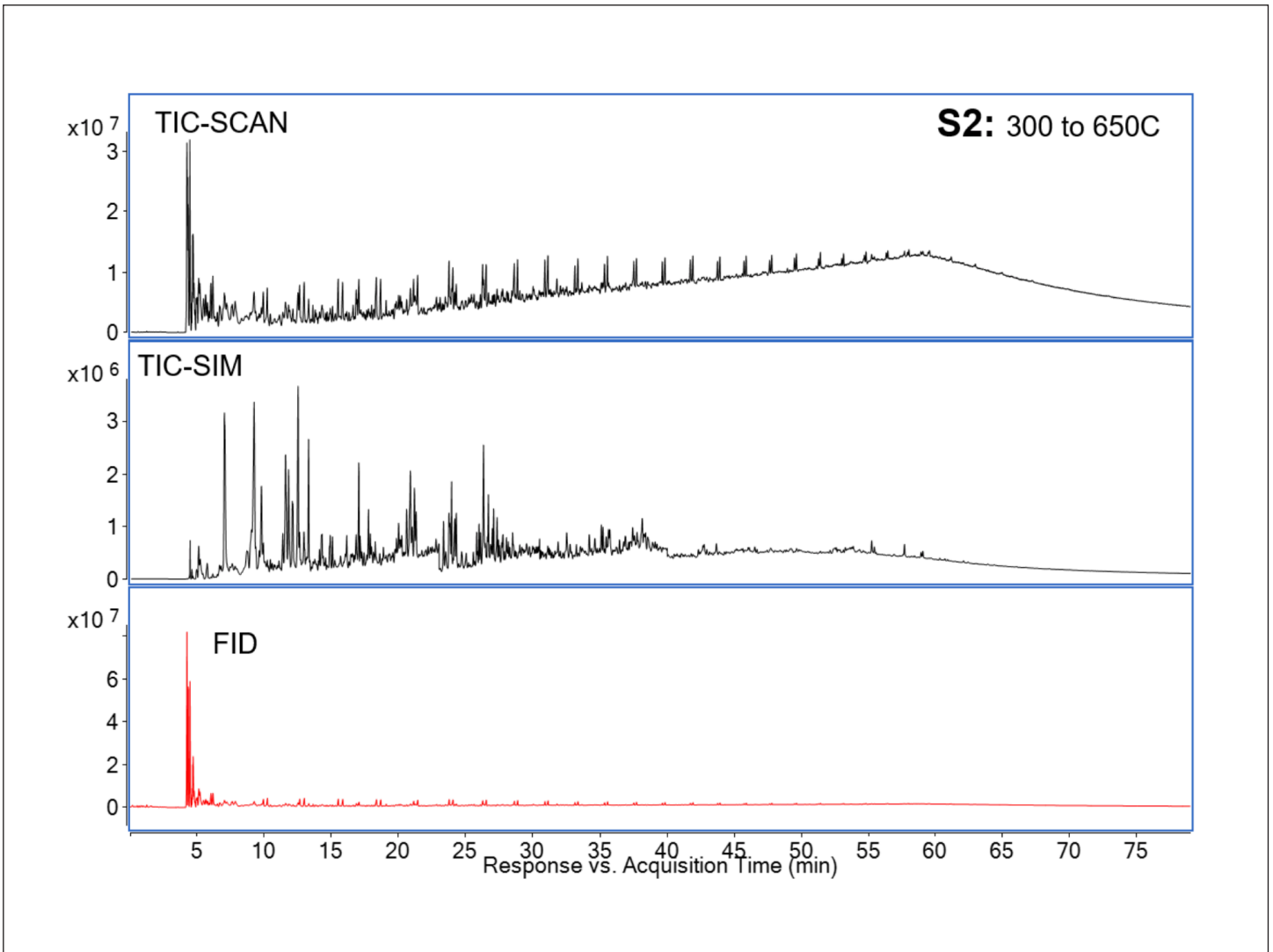


Figure 8. TIC-SCAN, TIC-SIM, and FID of S2.

Figure 9 shows the FID and the $m/z = 57$ (alkane) and $m/z = 55$ (alkene) peaks. The same fragmentograms for aromatics and biomarkers plotted for S1 may be

plotted for S2. In general, the distributions are quite different, especially for the biomarkers.

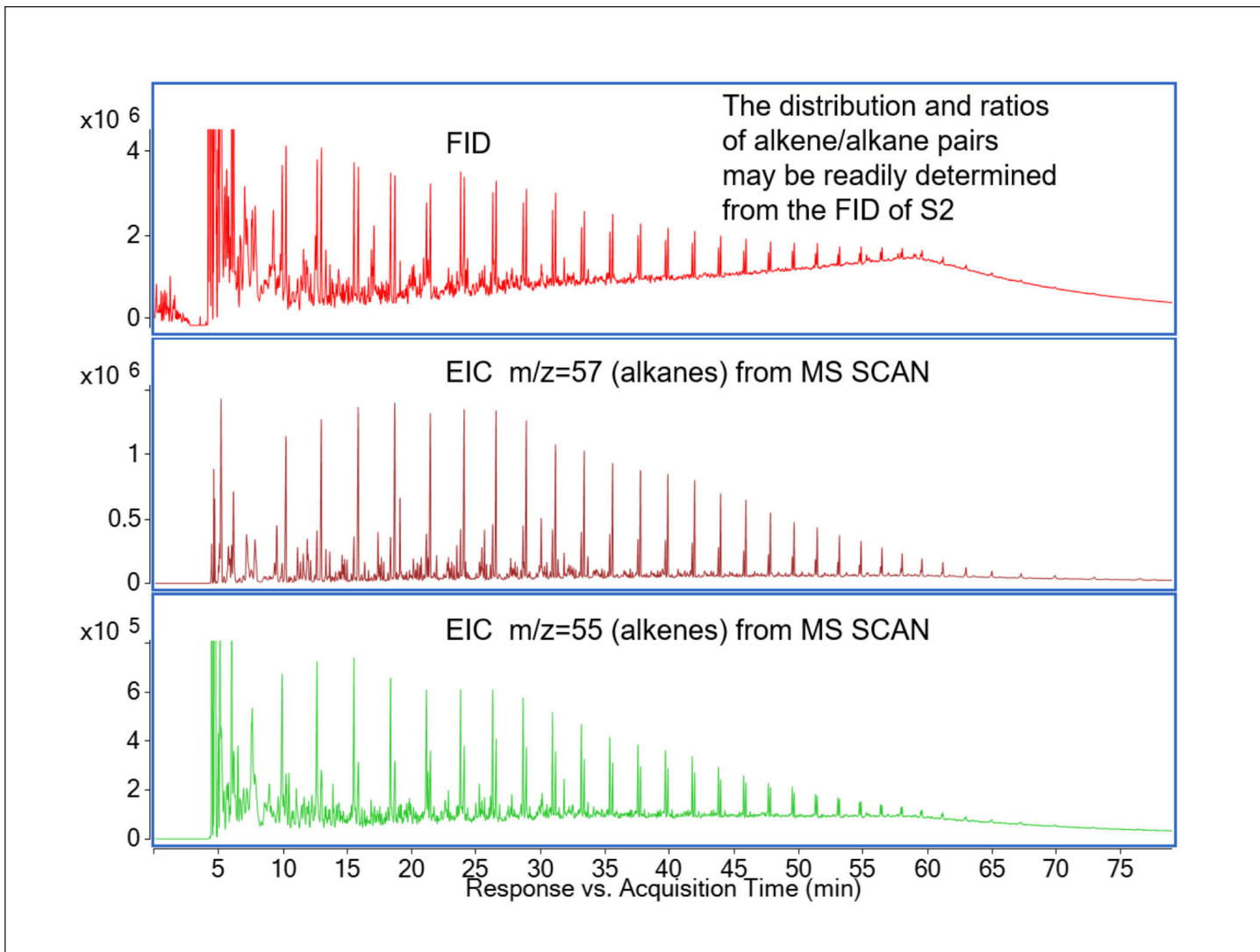


Figure 9. FID and EIC ($m/z:57$ and $m/z:55$) of S2.

Summary

In summary, the Thermal Vaporization/Pyrolysis-GC/MS technique using the vertical micro-furnace multi-mode pyrolyzer enables the analysis of S1, S2, and biomarkers present in a rock sample without any solvent extraction or sample preparation. The obtained detailed information from this technique can reveal critical facts about the recoverability and the potential for oil production.

Thermal extraction-GC/MS offers the geochemist a means of:

- ✓ Simplifying sample preparation.
- ✓ Improved accuracy of organic determination.
- ✓ Precise information on chemical species present.
- ✓ Ability to perform Rock Eval pyrolysis and biomarkers speciation on the same system.

About Quantum Analytics

Quantum Analytics is the authorized distributor for Frontier Laboratories in the United States and Canada. Contact us today to discuss your application needs, or to request a quote for the Frontier Multi-Shot Pyrolyzer.

Visit www.LQA.com/Pyrolysis to learn more.

Appendix A

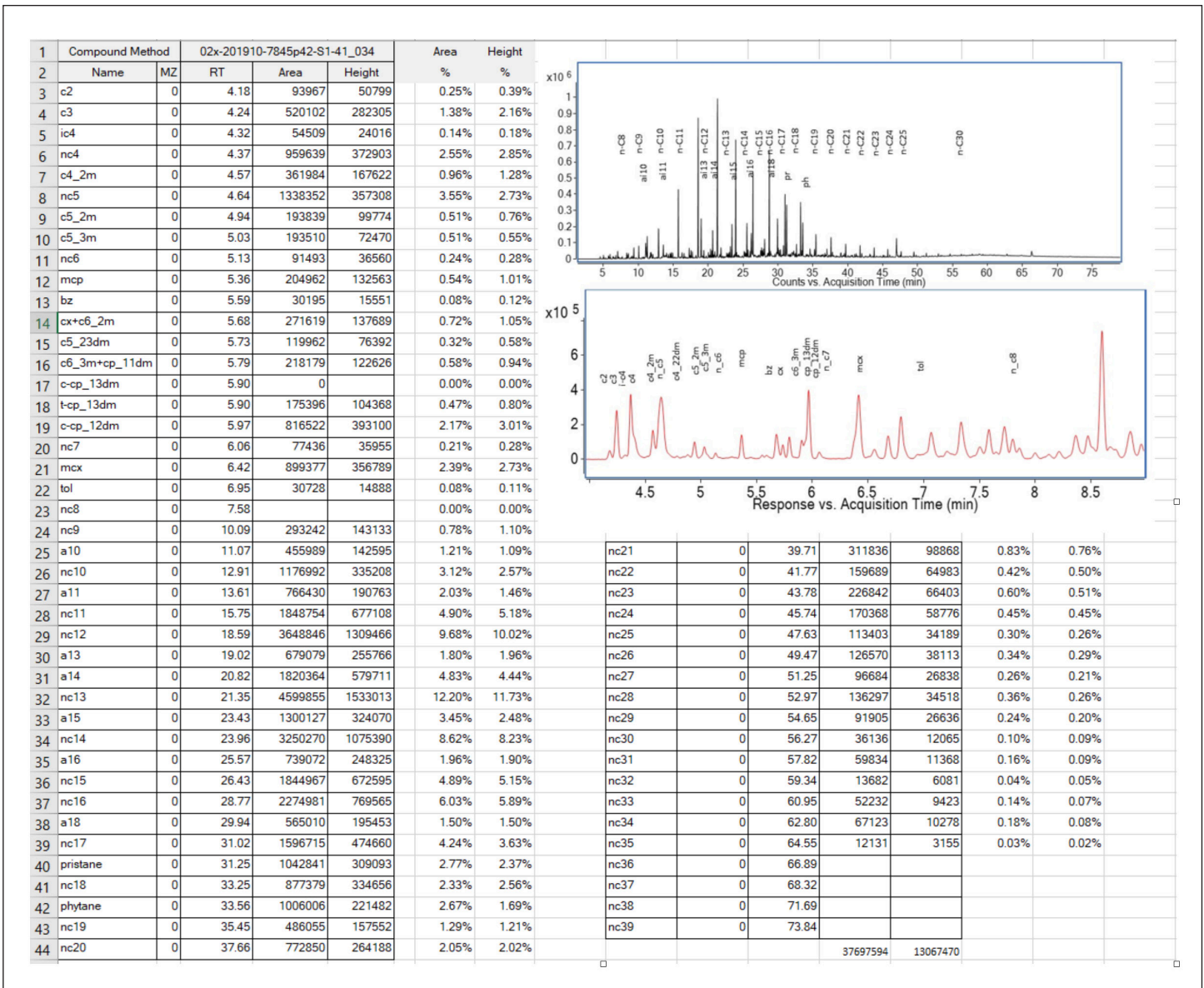


Table 1. Measurements of peak areas from the FID chromatogram, along with an expanded plot of the FID trace from 4 to 9 minutes.

Appendix B

Compound Method		201910-7845p42-S1-41			Area	Height		
Name	MZ	RT	Area	Height	Ratio	Ratio		
T19-C19 tricyclic terpene	191.2	36.14			Area Ratio=Area*100/Area of 30-Hopane			
T20-C20 Tricyclic terpene	191.2	38.00			Height Ratio=Area*100/Height of 30-Hopane			
T21-C21 tricyclic terpene	191.2	40.44	55646	17353	11.01	12.50		
T22-C22-tricyclic terpene	191.2	42.16	71042	20691	14.06	14.90		
T23-C23-tricyclic terpene	191.2	44.10	104704	30870	20.73	22.23		
T24-C24-tricyclic terpene	191.2	45.19	86277	27392	17.08	19.73		
T25S	191.2	47.33	42180	15049	8.35	10.84		
T25R	191.2	47.39	54777	15002	10.84	10.80		
TET24	191.2	48.78	20349	6194	4.03	4.46		
T26S-C26 Tricyclic Terpene(22S)	191.2	48.89	29910	9220	5.92	6.64		
T26R-C26 Tricyclic Terpene(22R)	191.2	49.03	31700	9512	6.27	6.85		
25TNH	191.2	54.57			0.00	0.00		
Ts-C27 17α, 21β-22, 29, 30-Trisnorhopane	191.2	54.57	58347	16815	11.55	12.11		
Tm-C27 17α, 21β-22, 29, 30-Trisnorhopane	191.2	55.22	79541	22977	15.74	16.55		
H28-C28 17α, 21β-28, 30-Bisnorhopane	191.2	56.73	29592	6810	5.86	4.90		
T28R-C28 Tricyclic-Terpene (22R)	191.2	52.64	36979	10855	7.32	7.82		
T28S-C28 Tricyclic-Terpene (22S)	191.2	52.40	33360	10936	6.60	7.88		
H29-C29 17α, 21β-Norhopane	191.2	57.45	251989	72290	49.88	52.06		
Ts29-C29 18α, 30-Norbornohopane	191.2	57.55	74250	23406	14.70	16.86		
NM-C29 17β, 21α-30-Norhopane (Normoretane)	191.2	58.19	26191	6649	5.18	4.79		
T29R-C29 Tricyclic-Terpene (22R)	191.2	53.71	46498	14162	9.20	10.20		
T29S-C29 Tricyclic-Terpene (22S)	191.2	53.42	45331	14079	8.97	10.14		
DH30-C30 Diahopane	191.2	57.79	3623	1551	0.72	1.12		
Ol-C30 18α-Oleanane	191.2				0.00	0.00		
H30-C30 Hopane	191.2	58.82	505202	138856	100.00	100.00		
M-C30 17β, 21α-Hopane (Moretane)	191.2	59.40	42933	9521	8.50	6.86		
G-C30 Gammacerane	191.2	61.17	69250	14523	13.71	10.46		
T30R-C30 Tricyclic-Terpene (22R)	191.2	55.64	47303	11474	9.36	8.26		
T30S-C30 Tricyclic-Terpene (22S)	191.2	55.32	41801	12188	8.27	8.78		
H31S-C31 17α, 21β-Homohopane (22S)	191.2	60.43	185426	47261	36.70	34.04		
H31R-C31 17α, 21β-Homohopane (22R)	191.2	60.62	128740	32913	25.48	23.70		
H32S-C32 17α, 21β-Bishomohopane (22S)	191.2	61.79	113442	26413	22.45	19.02		
H32R-C32 17α, 21β-Bishomohopane (22R)	191.2	62.10	78268	18475	15.49	13.31		
H33S-C33 17α, 21β-Trishomohopane (22S)	191.2	63.55	67936	15038	13.45	10.83		
H33R-C33 17α, 21β-Trishomohopane (22R)	191.2	64.03	46776	9846	9.26	7.09		
H34S-C34 17α, 21β-Tetrahomohopane (22S)	191.2	65.62	40841	7897	8.08	5.69		
H34R-C34 17α, 21β-Tetrahomohopane (22R)	191.2	66.29	24756	12806	4.90	9.22		
H35S-C35 17α, 21β-Pentahomohopane (22S)	191.2	68.01	30719	4880	6.08	3.51		
H35R-C35 17α, 21β-Pentahomohopane (22R)	191.2	68.95	13887	2516	2.75	1.81		

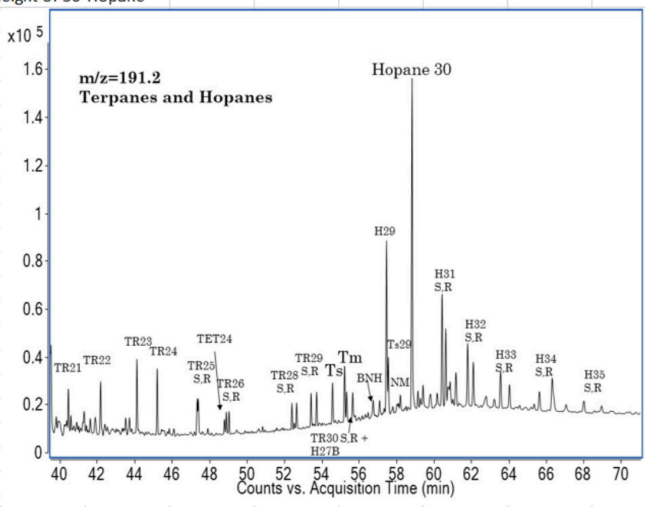


Table 2. Calculated results for the terpanes and hopanes relative to 30-Hopane normalized to 100.



About Quantum Analytics

Quantum Analytics is a value-added distributor of analytical equipment, offering of customized technical services, including cross-platform system integration, installation, training and support. Whether you are starting up a new lab or expanding your testing capabilities, we are here to help you take your science to new levels.



Instrumental to your success

 www.LQA.com |  LQA@LQA.com